


LabChart Overview

Data Acquisition and Analysis Software

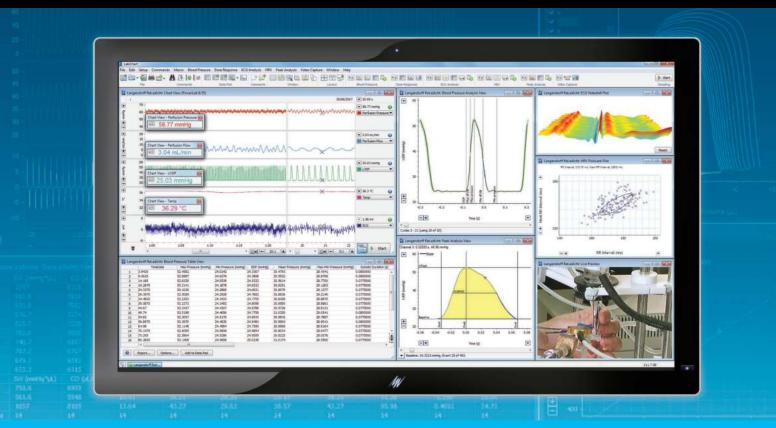

Designed specifically for life science, LabChart provides up to 32 channels for data display and analysis options that are powerful and easy to use.


What is LabChart?

LabChart data acquisition and analysis software creates a platform for all of your recording devices to work together, allowing you to acquire biological signals from multiple sources simultaneously and apply advanced calculations and plots as your experiment unfolds.

Quick to set up and simple to use, LabChart tracks every action you take and never modifies your raw data, ensuring the integrity of your results so you can focus on the true insights of your research.

Use LabChart to:

- Record data via PowerLabs and other compatible data acquisition devices
- Preview and optimize
 recordings
- Import and export data easily
- Annotate data with comments
- Automate recording and analysis tasks
- Convert raw data into useful units
- Change recording settings in seconds
- Recall data and experimental settings
- Generate customized stimulus outputs

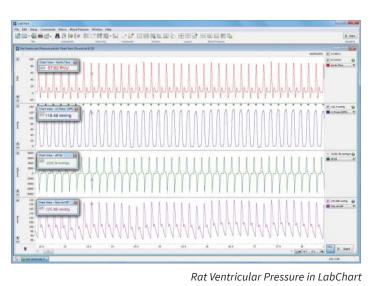


Streamlined Workflow

"Of most importance was the fact that all data was recorded on a timealigned beat-by-beat basis - this is vital for the calculation of transfer function analysis based on the relationship between cerebral blood flow and blood pressure."

Professor Ainslie Canada Research Chair Cerebrovascular Physiology



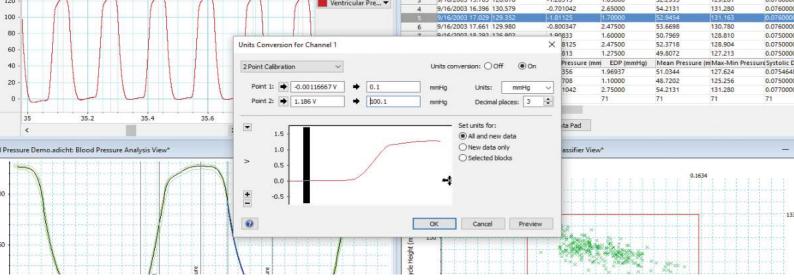

Signals

LabChart is flexible to the needs of novel research, with the ability to detect, display, and analyze a broad range of signals across many life science applications.

Supported signals include:

| Angles | EOG | Intracellular Recordings | Respiratory Gas Analysis |
|-------------------------------|--------------------------|--------------------------|--------------------------|
| Dissolved Gasses | Extracellular Recordings | NIBP | Sounds |
| dP/dT | Fluid Flow | рН | SPO ₂ |
| ECG | Force | Pressure | Temperature |
| EEG | Glucose | Pulse | Tissue Perfusion |
| Electrical Stimulation | GSR | RER | Video |
| EMG | Heart Rate | Respiratory Flow | Volume |

Research Applications

Designed specifically for life science researchers, LabChart provides all the calculations, reports, and plots you will need for human, animal and in vitro applications.

Animal Human Autonomic Autonomic Behavior, Sleep Cardiovascular and Neuroscience Exercise and Sport Cardiovascular Anima Neurophysiology Autonomi Telemetry Human Psychophysiology **Tissue and Circulation** Respiratory Sleep In Vitro In Vitro Speech Pathology Electrophysiology Tissue and Circulation **Isolated** Organ

Innovate and extend your studies

Providing the flexibility to extend your studies into new territories, it is easy to add new signals and measurements in LabChart as your research grows and takes on new directions. Looking to stand out from the crowd? LabChart also allows you to create your own advanced calculations and macros to customize your analysis.


If you have a particular research need you would like to discuss, get in touch with our experienced support team. We can work with you to customize an effective solution to record and analyze the specific data that you need.

LabChart Features

LabChart is designed for quick and easy use, allowing full control of your data. Advanced features stay out of the way until you need them, letting you do the basics with minimal fuss.



LabChart combined with A-M Systems hardware, displaying Spike Histogram analysis.

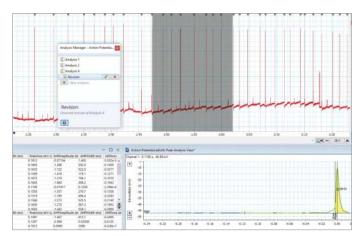


Analysis Manager

Save and label your selections and analysis regions, then recall them with a click when needed again.



Arithmetic

Build your own customized formulae, detection, and analysis within LabChart.

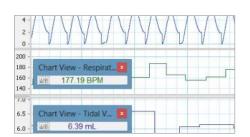


Cyclic Measurements

Combines flexible and clever cycle detection to create a simple to use analysis for periodic waveforms.

| Devices and Dravels | Besum and Chanas | | | a 🚺 | | |
|---|--|--------------------|---|--|--|--|
| Pudi Difer Deuton | Schate Ferendecte | | Name of damage | H (\$) | the ri | |
| Marker Rescaled Connection Marker Rescaled Connect Co | Denne 22 Proceeding ACD 35 22 Proceeding 207 9 | Ownsh 14 8-0 | Otennel R Dannel J R Dannel J R Dannel J R Dannel S R Dannel | Input Input 1 (Prove) Input 2 (Prove) Input 2 (Prove) Input 5 (Prove) Input 5 (Prove) Input 5 (Prove) Input 5 (Prove) Input 1 (Prove) Input 2 (Prove) Input 2 (Prove) Input 2 (Prove) | 4433 * 4433 * 4434 * 44344 * 44344 * 4434 * 4434 * 4434 * 4434 * 4434 * 4434 * | |
| Feature Manager Frequently Advections | 4 Although the | | | OK. | (ma | |

Device and Channel Discovery

Easily manage which devices, inputs and LabChart channels will be acquiring data - all from one panel.

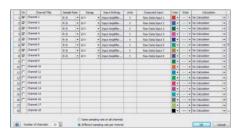


Channel Settings

Get an overview of, and easily change your recording settings, calculations and channels.

| AL S | 2.+ 000 | te]51 to te]513 | | | | WICH HAR I HAVE I |
|--------------|---|---|--|-------------------------------------|--|--|
| A | 6 Charmel 1 Avg Cvc Troquency Har | C Charriel 1 Average Cyclic Rate IRPM | D Channel 1 Maximum Value HV | E Charves 1 Taxe at Maxmum | F Channel 1 Maxmum Power V ^R | Ghasman I Maamum Power Frequency Hz |
| 1 0.01-51.0 | 234.984E | 14099.072 | 131,25 | 0:01:51.1351 | 0.0 | 175.7813 |
| 2 | Channel 1 | Channel 1 | Chantel 1 | Channel 1 | Chernel 1 | Channel 1 |
| 3 Time | Avg Freq | Avg Rate | Maximum | Max Time | Max Power | Max Power Frequency |
| 4 | Hz | BPM | mV. | | Va | Hz |
| 5 0 00 00 0 | 302.7852 | 18167 1132 | 395,7813 | 0.00.00.03238 | 0.0002 | 312.5 |
| 6 0.00.01.0 | 254.867 | 15292 022 | 321 1963 | 0.00.01.1457 | 0.0002 | 312.5 |
| 3 0.00.02.0 | 281.7287 | 16003 7199 | 266.5625 | 0.00.02.2916 | 0.0002 | 312.5 |
| 8 0.00 03.0 | 277.3466 | 16640.8135 | 308.8125 | 0.00.03.2871 | 0.0002 | 312.5 |
| 9-0.00.04.0 | 272.2276 | 16333 6542 | 312.25 | 0.00.04.84415 | 0.0003 | 312.5 |
| 10 0 00 05.0 | 249-4001 | 14964 1036 | 276.1875 | 0.00.05.1304 | 0.0002 | 312.5 |
| 11 0.00 06.0 | 274.2485 | 16454 9097 | 339.8125 | 0.00 06.7881 | 0.0002 | 312.5 |
| 12 0 00 07 0 | 292,5903 | 17555 0002 | 307.0625 | 0.00.07.9068 | 0.0002 | 312.5 |
| 13 0 00 08.0 | 288.9575 | 17237.4511 | 275 375 | 0.00.08.0461 | 0.0002 | 175,7813 |
| ** **** | ACC ADD. | ARTIGE MODE | And some | is one one remainst | 0.0000 | 242.6 |

Data Pad

Convert your time based data and calculations from chart view into discrete values for statistical analysis.

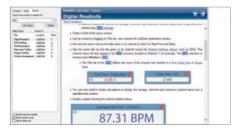


DVM

Display raw data or calculation signals as a large numerical display.



Comments

Seamlessly navigate your data using the annotations feature, either during or post data acquisition.

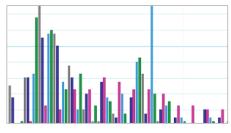


Data Plots

Visualize analysis easily with four different plot types and curve fits (Windows only).

| | 10 | | | - |
|--|-----|--------------------|----------|---|
| General | On | Product + | Download | |
| A | 12 | P/Loop New) | | |
| Installed Recommended Updates | 100 | Autio Dutout | | |
| Categories | 10 | Blood Pressure | | |
| Cardo-respiratory | 105 | Cardiac Output | | |
| Chemistry | 10 | Circadian Analysis | | |
| Data Visualization Bectrophysiology | 15 | DMT Normalization | | |
| File Translation | 83 | Dose Response | | |
| LabOhat Pro Modules | 5 | ECG Analysis | | |
| Measurement and Calibration New | 147 | Event Manager | | |
| Lainy | 10 | Export WAV | | |
| | 10 | HRV | | |


Feature Manager

Discover and install new features and updates, or toggle off features and menus you aren't planning to use right away.

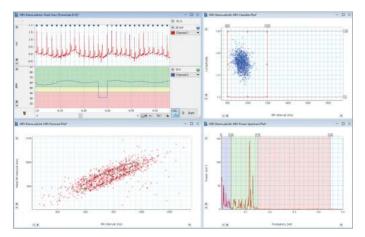

Guide Lines

Guide Lines provide a quick but powerful way to divide your data amplitude into regions of interest.

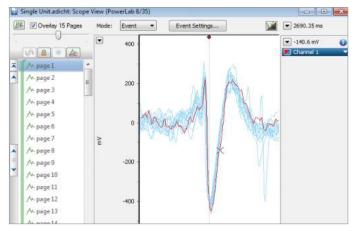

Scope View

Display, overlay, average and analyze periodic or evoked waveforms in real time.




Macros

Record your actions for repetitive operations, or get under the hood and write custom processes in VB Script.




Stimulator

Create a custom protocol to stimulate your preparation directly or to control a stimulus isolator.

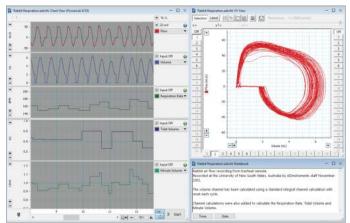

Layout

Display your screen the way you want. Smart Tile quickly snaps windows into a neat arrangement, or save your own layout preference.

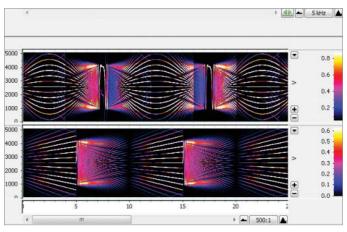

Spectrum

ile to play ba

O Loop file forever Reset file to beginning

8

device (Windows only).

Playback File

2 Point Calibration

150

50 0 -50 -100 -150

oint 1: 📦 -0.000378989

Units Conversion

Point 2: 🔿 6.60399 V

Playback File - ECG and Pulse.a | Browse...
Use sample rate from file
Include comments
Playback mode:
Stop at end of each block
Stop at end of file

OK Cancel

Stream pre-recorded data as if the data

Units conversion: 🗇 Off 🛛 🙆 On

Decimal places: 2

Set units for:

New data only

wHa

mmHg 🔹

were coming live from an attached


+ 100

This simple two point calibration tool

other common units of measurement.


makes it easy to convert volts into

Convert signals into component frequencies that can be extracted directly into data pad, or viewed as a power spectrum density, spectrogram or channel calculation.

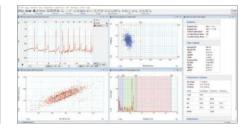

Split Screen

Show traces with two separate x-axis displays, ideal for viewing long term trends alongside latest data.

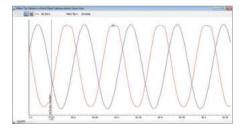


Zoom View

Take a closer look at an event, compare the timing between two overlaid channels, and export your images to publish.


LabChart Extensions

Extensions are optional features available with any version of LabChart and allow additional data formatting, filtering, visualization, measurement and calculation.

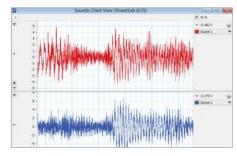


Audio Output

Allows data to be played back as sound. Useful for playing back input such as heart, animal and voice sounds.

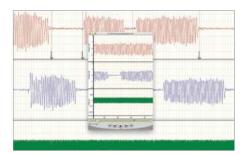


Export QuickTime

Save LabChart data files as QuickTime movies, then view on any supported platform or embed in other media.

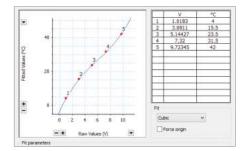


Multipoint Calibration Convert raw voltage signals from PowerLab into your desired units.

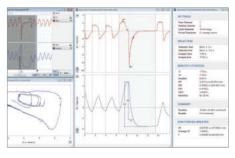


Spirometry

Derive respiratory parameters based on flow and volume such as VE, VT, PIF, PEF, FVC and FEV₁.

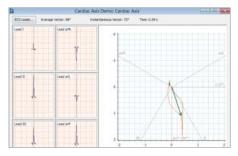

Cardiac Axis

Automate calculations of frontal plane ECGs and animate cardiac vector display.

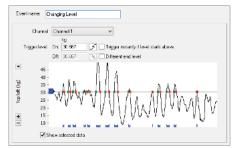


Event Manager

Monitors incoming signals, detects userdefined events and performs user-defined actions.


| Generate output when all | w of \$ | e following o | conditions are met; |
|-----------------------------|---------|---------------|---------------------------|
| Non 39071 - a lab | ₩£ + | 100000 | U D Holman I S |
| Adput settings | | | |
| III Paed width plates | 100.0 | 14 | must after pater starts 0 |
| Doley julin for | 0 | 1. | |
| Content washing of publics: | ÷ | | |

Fast Response Output

Control digital and analog outputs of supported PowerLab units.

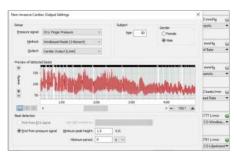

Non-Invasive Cardiac Output Allows a real-time estimation of continuous cardiac output to be calculated from an NIBP trace.

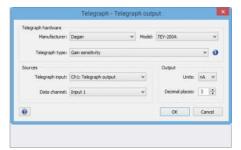

Telegraph

Make use of the gain telegraph output from an electrophysiological amplifier.

LabChart Remote App

Connect your iOS device to your sampling computer and operate certain functions of LabChart remotely via WiFi.

| Scheduled d | Sevices. | | Options | | | |
|-------------|-------------------|---------|---------|-------------|----------------|------|
| PowerLab | | 9 | | Autosave | Add comment | 8 |
| Recording a | etup | | | | | |
| Record: | 5.0 | ninutes | | Prom: | 09:00 a.n. 🚖 1 | oday |
| Every: | 60.0 | minutes | v | to: | 05:00 p.m. 🕀 | |
| | | | | () For: | 4 2 days | |
| | | | | C Every day | | |
| LebChart de | eta files | | | | | |
| Sounds | | | | | | |
| Vaual Ev | oked Potenitals | | | | | |
| Reaction | Time to Vioual Cu | | | | | |
| EDR Stw | rtle | | | | | |
| THRU | | | | | | |

Scheduler


Schedule recordings across one or more LabChart data files.



LabChart makes it easier to go wherever your research leads


Specialized analysis with LabChart Modules

Modules provide highly specialized data acquisition and analysis features for specific research applications. The full suite of modules are included with a LabChart Pro License, or can be purchased individually as needed. Some modules are not available on Mac.

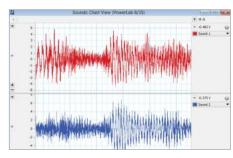


Dose Response

Generates dose response curves, EC50 values and additional parameters.

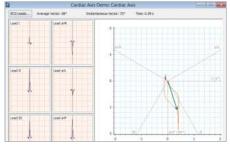


Metabolic

Provides real-time measurements of parameters such as VCO₂, VO₂, VE and RER.

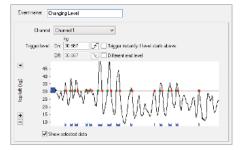


Blood Pressure

Automatically detects, analyzes and reports on parameters from arterial or ventricular pressure recordings.

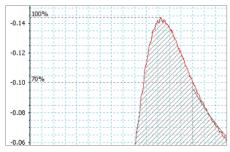


Spike Histogram

Detects, discriminates and analyzes extracellular spike activity generating a range of plots and statistics.

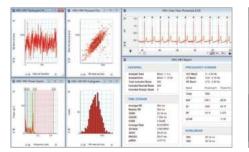


Heart Rate Variability

Displays and analyzes variation in the interval between heartbeats in human and animal ECG.

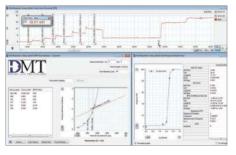


DMT Normalization

Calculates and standardizes optimal vessel pretension conditions using the wire myograph.

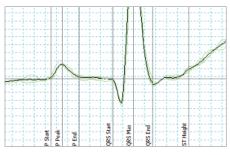


ECG Analysis

Detects and reports the onset, amplitude and interval times of PQRST from human and animal ECG signals.

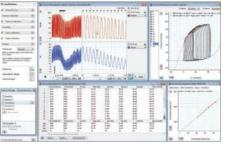

PV Loop

Analyzes left and right ventricular pressure/ volume data, calculates loop area plus a wide range of hemodynamic parameters.

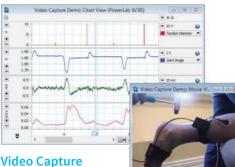


Video Capture Allows the synchronized

recording and playback of a Windows Media Video and LabChart data file.

Cardiac Output

Calculates cardiac output from a LabChart recording of a thermodilution curve measured in animals.

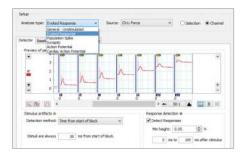

Peak Analysis

Automatic detection and analysis of multiple (but not overlapping) signal waveforms from recording.

Diverse Applications

"... the synchronized video is really a complete solution, taking swimming technique analysis to the next level."

Budd Termin Swimming Coach, TeamTermin



Hardware Compatibility

Signal conditioners and instruments

ADInstruments manufactures a wide range of software-controlled pre-amplifiers and offers auxillary devices for specialized investigations.

Transducers and accessories

• Built to last with 5 year warranty

We offer an extensive selection of instruments, transducers and accessories made by ADInstruments and world-leading brands.

• Simple to set up, USB connection to Windows or Mac

PowerLabC PowerLab

PowerLab is a high quality data acquisition device engineered to record precise, reliable, consistent data for research and education.

PowerLabs are capable of high speed sampling and form the core of many research data acquisition systems - they are compatible with instruments, signal conditioners, and transducers supplied by ADInstruments and many other leading brands.

- Automatically detected by LabChart
- High speed sampling (up to 200 kHz per channel)

PowerLab C + C Series Interfaces



PowerLab C is a digital data acquisition device that provides adaptive mains filtering, power management for peripheral devices (max 100W via USB-PD) and sub-µS time synchronization for up to four C Series compatible USB-C devices.

The C Series Front End Interface

converts analog data from ADInstruments Front-Ends such as Bridge Amps and Bio Amps to allow digital sampling by the PowerLab C. The **C Series Instrument Interface** provides 4 channels of input capability from any analog instrument to PowerLab C.

Research PowerLab Range

PowerLab 2/26 and 4/26

Available in either 2 and 4 channels, maximum sampling rate of 100kHz per channel, independent ADCs for each channel to keep data perfectly in sync.

LabChart Compatible Systems

Education PowerLab Range

PowerLab 15T and 26T Economically priced, high quality DAQ unit with built-in, fully isolated, twochannel bio amplifier, isolated stimulator, separate non-isolated analog inputs, and a differential-mode analog output.

As well as sampling from a broad range of instruments, signal conditioners and transducers via PowerLab, LabChart can also stream data directly from a range of compatible digital and wireless devices. These are available from manufacturers such as DSI, Oxford, DMT, Equivitial, Delsys and Kent.

LabChart for Education

Research-grade software for recording and analyzing life science data.

LabChart data acquisition and analysis software lets students gather biological signals from multiple sources simultaneously and apply calculations and plots as their experiments unfold.

Access a Range of Teaching Resources

To save you valuable time in creating course materials, we have developed over 150 LabChart exercises for education. Any of these exercises can be customized quickly and set up easily for your laboratory class.

Each lesson includes: Instructor Guides • Settings Files • Student Protocols • Preconfigured Lab Reports **Courses Covered:** Human Physiology • Animal Physiology • Psychophysiology • General Biology and Biochemistry Exercise Physiology • Neuroscience

Versions and Licenses

- Data Recording and Analysis

LabChart Pro

Recommended for researchers and educators who want the full range of modules and capability, including automation of specific measurements.

LabChart

Recommended for researchers and educators who want the capabilities without modules and automation of specific measurements.

Analysis only and collaboration


LabChart Pro Student License

Limited 6 month licence for undergraduate students. Provides full analysis, save and export functionality (no sampling). Ideal for students conducting data analysis of existing recordings within a curriculum.

LabChart Reader

Recommended for anyone who wants to view LabChart data on Windows or Mac. LabChart Reader doesn't allow saving of data - export image only. Has data pad for analysis.

| Main Features | LabChart Pro | LabChart | LabChart Pro Student License | LabChart Reader (Free) |
|---|--------------|--------------------------|---------------------------------|---------------------------|
| 5 years of updates + upgrades to all LabChart software | \checkmark | × | × | × |
| Open and read LabChart files | \checkmark | \checkmark | \checkmark | \checkmark |
| Analyze | \checkmark | \checkmark | \checkmark | \checkmark |
| Save and export | \checkmark | \checkmark | \checkmark | × |
| Record data | \checkmark | \checkmark | × | × |
| LabChart Modules | \checkmark | 🗴 (Purchased separately) | \checkmark | × |
| LabChart Extensions | \checkmark | \checkmark | \checkmark | × |

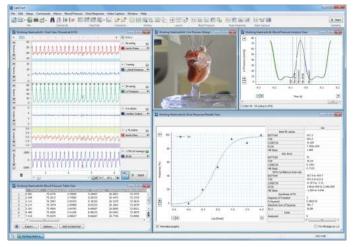

Radnoti Cardiovascular Blood Pressure analysis in LabChart

Maximize your potential

ADInstruments global training and support systems mean that there is always help at hand to support you with your endeavours and help you learn 'best practices' to move your research forward faster.

Data Integrity

"I like the fact that you can store all your data electronically. There is no need to create an auditing trail, because this is automatic. You can't remove bits of information. What is really interesting is that you can really go into the detail of the actual blood pressure trace and perform complex analyses very quickly. You simply export data into a standard template in Excel and have all your answers in seconds. Very smart."

Dr. Kengatharan,

President & COO, Armetheon, Inc.



ADInstruments equipment is used in the top 100 institutions for Life Science worldwide and is cited in more than 30,000 peer-reviewed papers.

PowerLab and LabChart are trademarks of ADInstruments Pty Ltd. All other trademarks are the property of their respective owners. Products supplied by ADInstruments are intended for use in research and teaching applications and environments only.

腾 <<p>(준 🚳 🛧 💽 😡 GLP

Visit adinstruments.com or contact your local ADInstruments representative for more information

Australia | Brazil | Europe | India | Japan | China | Middle East | New Zealand | North America | Pakistan | South America | South East Asia | United Kingdom



